

Laboratory Infrared Spectroscopy of atmospheric constituents

Background

Remote sensing for determination of the distribution of atmospheric trace gases relies on the spectroscopic line parameter database (e.g. HITRAN, SAO or JPL), such as line position, line strengths, pressure broadening or absorption cross sections. The database is mandatory to link the characteristic spectral signature obtained by remote sensing to the trace gas distribution in the atmosphere. Remote sensing of the atmosphere is carried out from ground-, airplane-, balloon- or satellite-based platforms. The spectroscopic database is usually obtained from laboratory experiments. The following sections introduce the infrastructure related to lab spectroscopy of DLR's Remote Sensing Technology Institute.

High Resolution FT Spectrometer

The Fourier Transform (FT) spectrometer at DLR is the commercial high resolution BRUKER IFS 125HR instrument covering the spectral range from 10 to 50000 cm^{-1} with a maximum resolution of 0.001 cm^{-1} (30 MHz). The instrument can be evacuated to avoid absorption of radiation by water and carbon dioxide in the ambient air. The spectrometer is equipped with a number of highly sophisticated absorption cells covering absorption paths from 0.15 m to 0.5 m and temperatures from 200 K to 1000 K. Several sensitive detectors exist covering the entire electromagnetic spectrum from FIR to UV. For the MIR region the detector optics was optimized to yield up to 10 times improved sensitivity over standard commercial detectors.

Multireflection cell

The multireflection cell is amongst others used for measurements of unstable atmospheric species which are produced in a steady state flow. Since only low average number densities are obtained, a large absorption path is required in order to provide a reasonable absorption for

the features of interest. Furthermore, column densities (number density times absorption path length) in the atmosphere, especially for stable species, e.g. water, can be very large, requiring also large column densities for the laboratory measurements. The large absorption path is realized by a setup applying spherical mirrors allowing the radiation to pass through the cell several times. The optical setup of the cell mirrors as well as the transfer optics for in- and out-coupling of the radiation from the FT spectrometer was designed by means of a special ray tracing program written at DLR. The cell was optimized for use in the FIR spectral region (10 – 500 cm^{-1}). The maximum absorption path length is 100 m, the base length of the cell 0.8 m and the diameter 0.2 m. Temperatures down to 200 K are achieved by circulating coolant through the double-jacket wall. The cell is unique with respect to its sensitivity for measurement of unstable species in the FIR. Therefore, FIR-FT spectra of species like OH, HO₂, BrO, ClO and Cl₂O₂ could be measured for the first time.

Coolable short path cell

The short path cell is used for measurements of stable atmospheric species at stratospheric temperatures down to 195 K. The cell exhibits an excellent temperature homogeneity (0.5 K, verified by spectroscopic gas temperature measurements) which is crucial for precision measurements of line strengths and pressure broadening parameters. The absorption path is 0.25 m with an inner diameter of 0.04 m. The window material can be selected according to the required spectral range and chemical stability, e.g. polyethylene for the FIR, or AgCl for the MIR and aggressive gases. The cell body consists of a Duran glass double jacket allowing to flow coolant through the entire length of the cell wall. Windows are mounted with good thermal conductivity to the



Bruker IFS 125HR



look into the instrument

Deutsches Zentrum
für Luft- und Raumfahrt e.V.
German Aerospace Center

Remote Sensing Technology Institute
Oberpfaffenhofen
D-32234 Weßling

Dr. Manfred Birk
Phone: +49 8153 – 28 3084
Fax: +49 8153 -28 1337
Manfred.Birk@dlr.de
www.DLR.de/eoc



cell body to further improve temperature homogeneity. The cell has been applied for investigation of the molecules O_3 , N_2O_5 , $ClONO_2$ for atmospheric remote sensing.

Heatable short path cell

This cell is used for measurements under elevated temperatures up to 1000 K. The absorption path is 0.16 m with an inner diameter of 0.04 m. The cell body is made of quartz glass, the windows are CaF_2 , providing high temperature resistance combined with mechanical stability while at the same time allowing MIR measurements down to 1000 cm^{-1} . The cell is heated electrically by wires attached to the cell body and the gas inlet tube. Special care was taken to keep the windows at approximately the same temperature as the cell to improve temperature homogeneity. The cell is equipped with in- and outlet tubes to allow for a slow flow through the cell in case species with limited lifetime at high temperatures are measured. Line parameters or absorption cross sections of CO , CO_2 , H_2O , NO , NO_2 have been measured up to 1000 K applying this cell in support of non-intrusive measurements of aircraft exhaust.

Production of unstable atmospheric compounds

Several atmospheric trace gases are very instable, reactive (mainly radicals) and cannot be bottled nor trapped. Of course such gaseous compounds cannot be maintained in an absorption cell statically but they can be produced in a flow reactor in steady state and pumped through an absorption cell. Between in- and outlet a concentration gradient is achieved invariant in time due to decomposition. The average particle density is low and will decrease with decreasing lifetime. Thus, a multireflection cell with a long absorption path length is required for FT-measurements. In order to obtain line

strength data from measured spectra the average particle density is needed. In contrast to stable compounds the particle density cannot be deduced from simple pressure measurements. However, in case of pure rotational transitions located in the FIR the line strength can be calculated from the permanent electric dipole moment of the molecule which is precisely known from Stark measurements. Thus, from FIR measurements the particle density can be derived allowing to obtain line strength data in other spectral regions. Commonly, the precursors for radicals are atoms of elements like fluorine, chlorine, oxygen or hydrogen which are produced in a microwave discharge in our case. The atomic species are produced in a steady flow reactor with other precursors. The gas flows which are almost entirely inert gas are measured and regulated by means of electronic mass flow controllers. The minimal residence time in the absorption cell is 0.5 s, the total pressure ranges from 5 to 20 hPa. So far the radicals OH , HO_2 , BrO and ClO have been measured, all of these being of atmospheric relevance.

Precision intensity measurements

The extension of the spectroscopic database of atmospheric trace gases requires accurate measurement of spectral intensities and line shapes. Whereas the accurate determination of line positions from FT spectra can be achieved easily the measurement of intensities is hampered by numerous systematic error sources. These error sources have to be characterised and minimised. Unfortunately, a line intensity standard is not available for the MIR region. However, line strengths and thus line intensities can be calculated accurately in case of pure rotational transitions, which can be observed in the FIR region. Thus, the FT measurement of the pure rotational transitions of CO

allows accessing and quantifying the systematic error sources. For the first time, line strengths were experimentally determined with accuracy better than 1 %. Line shapes together with line shifts are temperature dependent and measurements with specified uncertainties are very rare. Measurements of O_3 in the MIR and H_2O in the MIR/NIR have been performed.

References

- Jacquinet-Husson, N.; Scott, N.A.; et al. (Birk, M.; Schreier, F.; Wagner, G.) (2008) The GEISA spectroscopic database: Current and future archive for Earth and planetary atmosphere studies. *J. Quant. Spectroscopy And Rad. Transfer*, 109 (6), 1043-1059
DOI: 10.1016/j.jqsrt.2007.12.015
- Coudert, L.H.; Wagner, G.; Birk, M.; Baranov, Y.I.; Lafferty, W.J.; Flaud, J.-M. (2008): The $H_2^{16}O$ molecule: line position and line intensity analyses up to the second triad. *J. Molec. Spectroscopy*, 1 (251), 339-357
- Wagner, G.; Birk, M.; Gamache, R.R.; Hartmann, J.-M. (2005): Collisional parameters of H_2O lines: effect of temperature. *J. Quant. Spectroscopy and Rad. Transfer*, 92, 211-230
DOI: 10.1016/j.jqsrt.2004.07.023
- Rothman, L.S.; Jacquemart, D. et al. (Birk, M.; Wagner, G.) (2005) The HITRAN 2004 molecular spectroscopic database. *J. Quant. Spectroscopy and Rad. Transfer*, 96, 139-204
DOI: 10.1016/j.jqsrt.2004.10.008
- Rothman, L.S.; Gordon I.E. et al. (Birk, M.) (2009) The HITRAN 2008 molecular spectroscopic database. *J. Quant. Spectroscopy and Rad. Transfer*, 110, 533-572
DOI: 10.1016/j.jqsrt.2009.02.013

Deutsches Zentrum
für Luft- und Raumfahrt e.V.
German Aerospace Center

Remote Sensing Technology Institute
Oberpfaffenhofen
D-32234 Weßling

Dr. Manfred Birk
Phone: +49 8153 – 28 3084
Fax: +49 8153 -28 1337
Manfred.Birk@dlr.de
www.DLR.de/eoc